


PROPER USE OF FUME HOODS




DO




Understand the chemicals in use and refer to SDS if uncertain

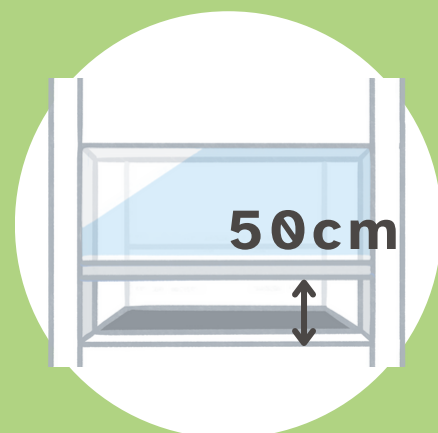

Ensure the sash is at 50 cm or lower

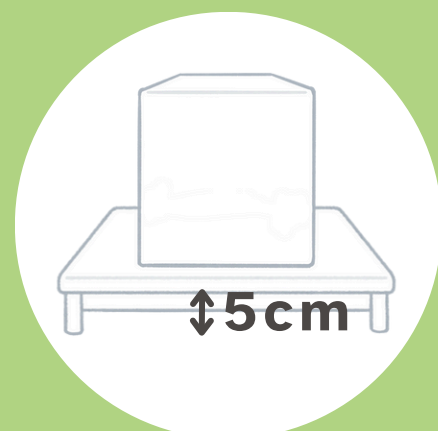

Raise large equipment at least 5 cm above the base to prevent airflow blockages

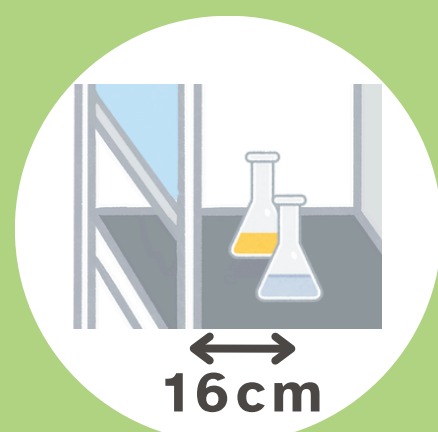

Keep all materials inside the hood at least 16 cm away from the sash opening



Ensure the hood is in operation and without flow alarm



Place the heat generating equipment near the back of the hood

DON'T



Put your head inside the fume hood



Store chemicals permanently in the fume hood



Leave the sash open when it is not in use



Use the fume hood for unintended purposes



Use any electronic extension cord inside the fume hood



Obstruct the obstacle detection sensor

